

# LAARMANN<sup>®</sup>

*Innovators in Solids*

## CEMENT CLINKER LM1000 PULVERISER RING MILL



Laarmann Group B.V.  
Op het Schoor 6  
6041 AV Roermond  
Netherlands

[info@laarmann.eu](mailto:info@laarmann.eu)  
[www.laarmann.eu](http://www.laarmann.eu)  
[webshop.laarmann.eu](http://webshop.laarmann.eu)  
+31 6 23 40 00 33

# CEMENT CLINKER APPLICATION

## Grinding parameters

Grinding system:	LM1000 Pulveriser Ring Mill
Accessories:	Grinding bowl 125ml / Tungsten carbide
Speed:	1400 rpm
Sample material:	Cement Clinker
Product feed size:	Pieces of <b>up to 16mm size</b>
Grinding time:	60 seconds
Sample amount:	51,49
<b>Result:</b>	97% < 63 micrometer after 60 seconds grinding time



## Recommendation

For the sample preparation of Cement clinker we recommend our Pulverizer Type LM1000 under the above mentioned conditions. A prolongation of the grinding time will improve the end-fineness. **The Liftboy** system for the LM1000 ensures an easy handling of the bowl without any manual action necessary.