



Cement clinker

Application

Market segment:	Construction Materials
Material:	Cement clinker
Feed size:	< 8 mm
Feed quantity:	80 g / sample
Material specification(s):	Dry, brittle
Customer requirement(s):	< 150 µm
Subsequent analysis:	X-ray Fluorescence Analysis

Solution

Selected instrument(s):	LM2000 Pulveriser Ring Mill
Configuration(s):	Grinding set hardened steel 100 ml
Working settings:	Revolution speed 1400 rpm
Time:	1 min.
Achieved result(s):	< 150 µm
Remark(s):	
Recommendation(s):	For this application the LM2000 Pulveriser Ring Mill is suitable under the above mentioned conditions.



Clinker

Application notes can only be used as a guideline. Each type material of has its own characteristics and therefore the outcome for each test may vary per material. We cannot accept any legal claims derived from this application note. This document is subjected to technical modification and errors.