You Tube Human Hair	
Application	
Market segment:	Biology
Material:	Hair washed and dried
Feed size:	Up to 40 mm
Feed quantity:	10 g
Material specification(s):	Fibrous
Customer requirment(s):	< 200 µm
Subsequent analysis:	Verification of drugs consumption / medical analyses
Solution	
Selected instrument(s):	LM PBM-1 Planetary Ball Mill
Configuration(s):	Grinding jar 250 ml stainless steel, 12 grinding balls Ø 20 mm stainless steel
Working settings:	Revolution speed 580 U/min.
Time:	60 min.
Achieved result(s):	< 200 µm
Remark(s):	
Recommendation(s):	For fine grinding of different hair samples the LM-PBM1 Planetary Ball Mill is suitable under the above mentioned conditions.

LAARMANN® **INNOVATORS IN SOLIDS**



Hair

Application notes can only be used as a guideline. Each type material of has its own characteristics and therefore the outcome for each test may vary per material. We cannot accept any legal claims derived from this application note. This document is subjected to technical modification and errors.