

LAARMANN[®]

Innovators in Solids

MAIZE CM1000



Laarmann Group B.V.
Op het Schoor 6
6041 AV Roermond
Netherlands

info@laarmann.eu
www.laarmann.eu
webshop.laarmann.eu
+31 6 23 40 00 33

MAIZE APPLICATION

Grinding parameters

Grinding system: Cutting Mill Type CM1000

Bottom sieve: 12,00 mm

Grinding time: 20 secs. each

Sample amount: 2,5 kgs

Result after sieving: Maize

100% < 6,3mm

65% 2,6 – 6,3mm

35% < 2,5 mm



Maize original sample



Amount app. 2,5 kgs



Picture 3 Processed sample

Recommendation

In order to grind maize samples we recommend the use of the CM1000 Cutting Mill under the mentioned conditions. Considering the above parameters the total capacity per hour is about 400 kgs per hour provided the feeding and sample extraction is continuously.