

LAARMANN[®]

Innovators in Solids

STRAW GRASS CM500 CUTTING MILL



Laarmann Group B.V.
Op het Schoor 6
6041 AV Roermond
Netherlands

info@laarmann.eu
www.laarmann.eu
webshop.laarmann.eu
+31 6 23 40 00 33

STRAW GRASS APPLICATION

Grinding parameters

Grinding system: Cutting Mill Type CM500

Accessories: Bottom sieves of 4mm and 1,5mm

Sample material: Straw

Product feed size: Pieces of **up to 40 mm length / Qty. 270 gram**

Grinding time: 1 minute

Result: Homogeneously milled



Original samples No 1. Medica



Processed Sample after CM500 grinding with 4,00 mm bottom sieve.



Processed Sample after CM500 grinding with 1,50 mm bottom sieve.



Original sample No 2. Straw



Processed Sample after CM500 grinding with 4,00 mm bottom sieve.



Processed Sample after CM500 grinding with 1,50 mm bottom sieve.

Recommendation

In order to grind and to homogeneously mix fibrous samples like Straw, grass etc. we recommend the use of the CM500 Cutting Mill under the above mentioned grinding conditions.