

LAARMANN[®]

Innovators in Solids

VARIOUS DRIED NUTS MICRO BALL MILL



Laarmann Group B.V.
Op het Schoor 6
6041 AV Roermond
Netherlands

info@laarmann.eu
www.laarmann.eu
webshop.laarmann.eu
+31 6 23 40 00 33

Product description:

Evaluation of particle size distribution



VARIOUS DRIED NUTS APPLICATION

- Grinding system: Lab Wizz 320 Micro Ball Mill
- Accessories: Stainless grinding jar 100ml/ grinding ball 30mm
- Sample material: Dried Kola nut, Tormentillae Rhizana, Cynara Scolymus
- Product feed size: Pieces of **up to 30mm**
- Grinding time: 30 seconds for sample 1.
60 seconds for sample 2 and 3

Recommendation

In order to grind the above mentioned sample materials we recommend the use of the Lab Wizz under the above mentioned grinding conditions.