


LAARMANN[®]

Innovators in Solids

MARBLE APPLICATION LMC200 JAW CRUSHER

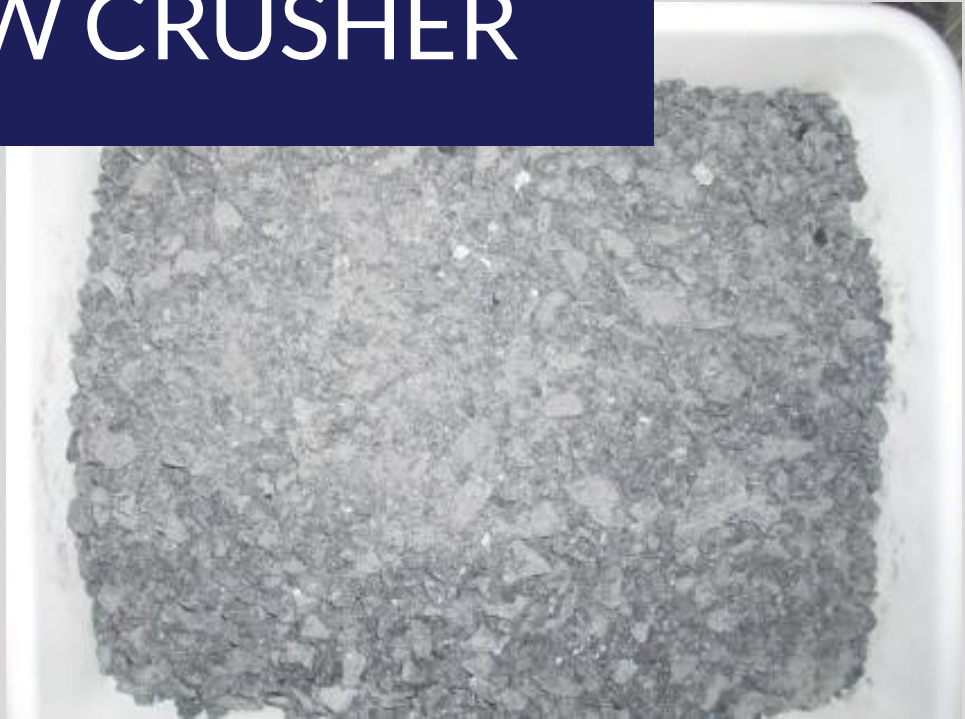

POLISILICON APPLICATION

Sample Material: Marble



Original sample marble



Final prepared sample

Grinding parameters:

Grinding system: LMC200 Jaw Crusher

Accessories: Breaking jaws and side liner of manganese steel

Sample material: Marble

Product feed size: 200mm length and 70mm diameter

Grinding time: 2 minutes per drill core.

Recommendation

For the sample preparation of a.m. material and the requested end fineness of 10-20mm our LMC200 Jaw Crusher is perfectly suitable but the material should be prepared in a wet condition to avoid agglomeration.